

Made
in
Germany



The **TDO Sahara** is a convection oven for fast and gentle deparaffining and drying of slides in the laboratory.

From the outside, the **TDO Sahara** offers full frontal view into the drying chamber through the generous glass window. This gives the user full control over the current state of the slide inside at all times. A magnetic locking mechanism ensures that the warm air remains inside the device. This way, the temperature remains stable and the energy expenditure is not unnecessarily increased.

The **TDO Sahara** warms up to operating temperature in just about a minute. The drying times are set using a touch control element. The user can choose between a chosen drying time or continuous operation, which is particularly useful when a large number of samples is processed. Every specified drying cycle is followed by a cooling phase. Routine slide drying usually only takes a few minutes.

The drying chamber offers space for up to 48 baskets with 30 microscope slides, distributed on two stainless steel grids. On the bottom of the drying chamber there is a drip tray for safely collecting dripping paraffin or other liquids. Both grids and the drip tray can be easily removed for cleaning.

In addition to the magnetic locking mechanism that prevents fumes from escaping to the outside, an activated carbon filter helps to reduce the pollution of the laboratory air by solvent vapors that can arise inside the drying chamber during operation.



Technical Data TDO Sahara		Cat. No.
Drying Time:	adjustable from 1 min. to 99 hrs. 59 min., or endless operation	
Temperature:	20 °C to 80 °C, adjustable in 1 °C steps	
Dimensions Drying Chamber (W/D/H)	414 x 273 x 280 mm	
Dimensions Overall (W/D/H):	445 x 368 x 480 mm	
Weight:	ca. 23 kg	
Power Supply:	230 V / 50 – 60 Hz / 1250 VA	02-6800-00
	115 V / 50 – 60 Hz / 1250 VA	02-6900-00

Accessories	Purchasing Unit	Cat. No
Activated Carbon Filter (disposable)	1 pc.	67-1006-00



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
Distributed by:

Axlab A/S

+45 35 43 18 81

axlab@axlab.dk

www.axlab.dk

AxlabYOUR VISION.
OUR INSPIRATION**MEDITE** **MEDITE Medical GmbH**

Wollenweberstr. 12 • 31303 Burgdorf • Germany

Phone +49 5136 88 84-0 • Fax +49 5136 88 84-55

info@medite.de • www.medite.de

Subject to change without prior notice. Pictures and colours may vary.